




# The Universal Goniometer: Complete Data Sets on All Crystals, No Compromise

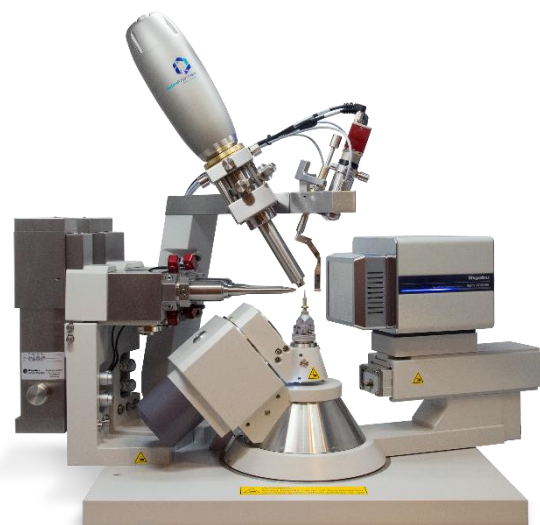

## Introduction

Based on the Universal Goniometer, the Rigaku Oxford Diffraction (ROD) XtaLAB Synergy X-ray diffractometers represent new and innovative tools for X-ray diffraction analysis in the home laboratory. The series include a single/dual wavelength microfocus sealed tube (the XtaLAB Synergy-S), and a microfocus rotating anode (the single wavelength XtaLAB Synergy-R and dual wavelength XtaLAB Synergy-DW). The central component for all ROD systems is the Universal Goniometer, a flexible and efficient full 4-circle kappa goniometer featuring the following:

- $\omega$  rotation speed of up to  $10^\circ/\text{sec}$
- $\omega$  rotation both in clockwise and counterclockwise directions to minimize goniometer repositioning time in between scans
- Minimum detector distance of 32 mm and maximum  $2\theta$  position of  $107^\circ$  to capture the highest resolution diffraction reflections
- Telescopic  $2\theta$  arm allowing for a maximum detector distance of 250 mm for large unit cell dimensions
- $2\theta$  and  $\kappa$  motion in both positive and negative directions and a  $\kappa$  range of  $\pm 179^\circ$  for the shortest data collection strategies

To demonstrate the flexibility and high performance of the Universal Goniometer, we present here the results from data collected on a small crystal of chlorothiazide crystallized in space group P1.

## Experimental Overview

A small crystal of chlorothiazide, with dimensions  $0.019 \times 0.035 \times 0.046$  mm, was captured within a thin layer of paratone oil inside a fiber loop mounted on a magnetic base (Figure 1). The crystal was mounted on the Universal Goniometer of a microfocus sealed tube XtaLAB Synergy-S X-ray diffractometer and data were collected at 100 K using ROD software, CrysAlis<sup>Pro</sup> 1.

**Table 1: XtaLAB Synergy-S specifications.**

<b>X-ray source</b>	PhotonJet-S Cu source with continuously variable divergence slit Beam FWHM = 110 $\mu\text{m}$
<b>Operating power</b>	50 kV x 1 mA = 50 W
<b>Goniometer Detector distance range</b>	4-circle Kappa with telescoping 2Theta arm 32 – 250 mm
<b>Detector Active area Frame rate Readout speed Pixel size Cooling</b>	Hybrid photon counting HyPix-6000HE 77.5 x 80.3 mm <sup>2</sup> Up to 100 Hz 0 ms in ZeroDeadTime mode 100 $\mu\text{m}$ air-cooled



**Figure 1: View of the chlorothiazide crystal mounted on the goniometer.**

## Data collection

A target completeness of 100% and redundancy of 3 at 0.837 Å on the Friedel mates merged was used to calculate the strategy in CrysAlis<sup>Pro</sup>. 26 scans were calculated, 9 at  $2\theta = -47.20^\circ$ , 4 at  $2\theta = -86.25^\circ$  and 17 at  $2\theta = -106.49^\circ$ . The exposure time was estimated by CrysAlis<sup>Pro</sup> for a target  $\langle I/\sigma(I) \rangle$  of 30 at 0.837 Å for the data unmerged, based on the  $\langle I/\sigma(I) \rangle$  calculated upon screening. The full data collection time was 2 hours and 51 minutes. Table 1 shows the data collection parameters.

Table 1: Data collection parameters.

<b>Generator settings</b>	50 W (50 kV, 1 mA)
<b>Wavelength</b>	1.54184 Å
<b>Temperature</b>	100 K
<b>Detector distance</b>	34 mm
<b>Low resolution exposure time</b>	1 seconds
<b>High resolution exposure time</b>	4 seconds
<b>Scan width</b>	0.5 °
<b>Total frames</b>	3176
<b>Total time</b>	2 hours 51 minutes

With exposure times of 1 second and 4 seconds per 0.5 ° for the low resolution and high resolution scans, respectively, diffraction from the crystal yielded small, very sharp reflections across the resolution range (Figure 2).

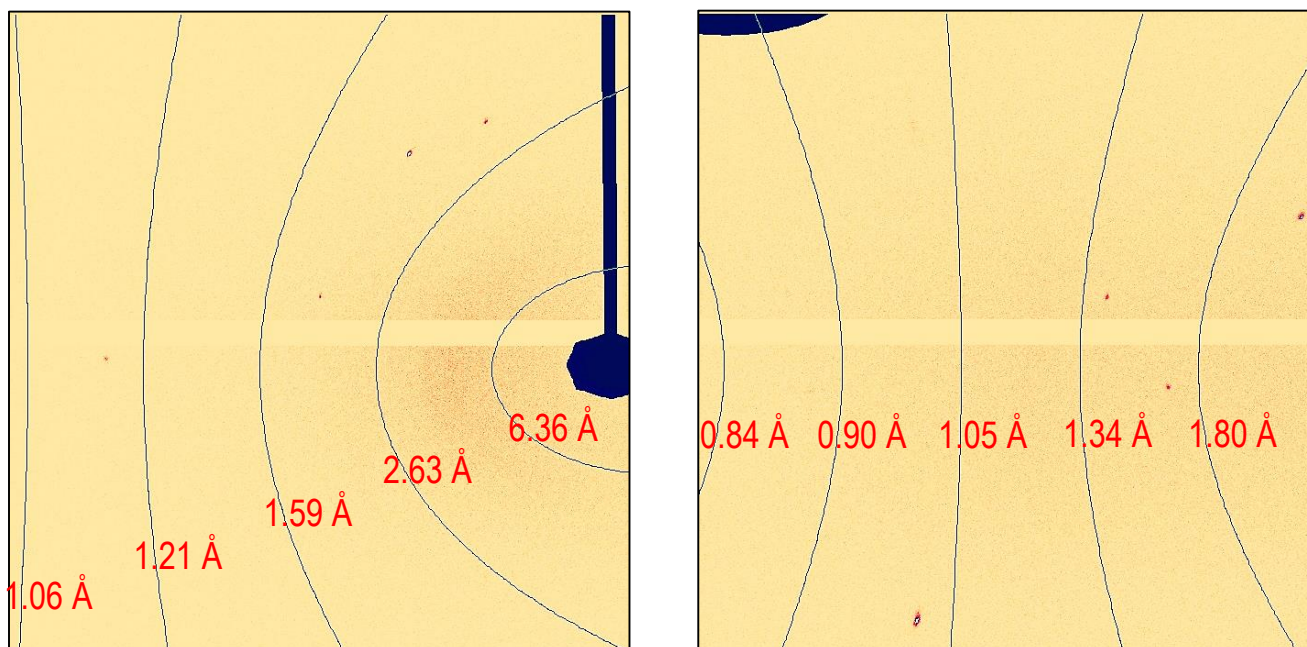


Figure 2: Example of diffraction images obtained at low resolution with 1 sec / 0.50° (left) and at high resolution with 4 sec / 0.50° (right).

## Results

Upon processing with CrysAlis<sup>Pro</sup>, the data set was 100% complete, including in the last resolution shell to 0.837 Å (Table 2). Data processing yielded an overall  $R_{\text{int}}$  of 3.6% and  $\langle I/\sigma(I) \rangle$  of 31.9 to 0.837 Å. The crystal structure (Figure 3) was automatically solved by ShelxT<sup>2</sup> in space group P1 and refined with ShelxL<sup>3</sup> within the AutoChem<sup>4</sup> interface. Hydrogen atoms were placed as riding atoms. Excellent structural statistics were obtained, with  $R_1$  at 2.51% and GooF at 1.070.

Table 2: Experiment and refinement details for the Friedel pairs unmerged.

Space group	P1
Unit cell	4.8469(3), 6.3804(3), 8.9211(4) Å 74.317 (4), 83.926(5), 80.475(5) °
Resolution	0.837 Å
Completeness (last shell)	100% (100%)
Redundancy (last shell)	3.1 (1.8)
$\langle I/\sigma(I) \rangle$ (last shell)	31.9 (20.1)
$R_{\text{int}}$ (last shell)	3.6% (5.8%)
Final R factors [ $I > 2\sigma(I)$ ]	$R_1 = 2.51\%$ $wR_2 = 6.29\%$
Goodness of fit	1.070
Largest residual peak/hole ( $\text{e}\text{\AA}^{-3}$ )	0.329 / -0.268

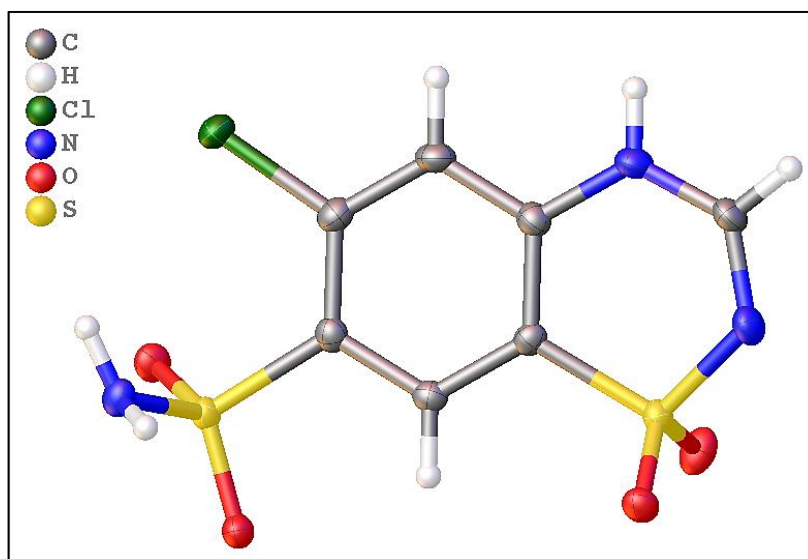


Figure 3: Representation of the model obtained for the chlorothiazide molecule after refinements to 0.837 Å.

## Conclusion

A 100% complete data set, both overall and in the last resolution shell to 0.837 Å, was collected on a small crystal of chlorothiazide (0.019 x 0.035 x 0.046 mm<sup>3</sup>) in under 3 hours, using the microfocus sealed tube X-ray generator XtaLAB Synergy-S configured with a 4-circles kappa Universal Goniometer. The crystal structure of chlorothiazide was solved readily and yielded excellent statistics. Such results demonstrate that:

- The Universal Goniometer allows for the collection of complete data sets, even in the lowest symmetry possible.
- Combined with the powerful strategy algorithm in CrysAlis<sup>Pro</sup>, the Universal Goniometer enables the collection of the best data sets in a minimum amount of time.

## References

1. Rigaku Oxford Diffraction, (2019), CrysAlis<sup>Pro</sup> Software system, version 1.171.40.45a, Rigaku Corporation, Wroclaw, Poland.
2. Rigaku Oxford Diffraction, (2019), AutoChem 4.0 software system in conjunction with Olex2 1.3, Rigaku Corporation, Wroclaw, Poland
3. Dolomanov, O.V., Bourhis, L.J., Gildea, R.J., Howard, J.A.K. & Puschmann, H., *J. Appl. Cryst.* (2009), **42**, 339-341.
4. Sheldrick, G. M. (2015), *Acta Cryst.* **A71**, 3-8.

## Rigaku Oxford Diffraction

9009 New Trails Drive, The Woodlands, TX 77381-5209

Tel.: (281) 362-2300 | FAX: (281) 364-3628 | [www.Rigaku.com](http://www.Rigaku.com) | [info@Rigaku.com](mailto:info@Rigaku.com)